

[View on rigaku.com](https://www.rigaku.com)

# EDXRF1631 - Zirconium on Aluminum



## Scope

This application note demonstrates the measurement of zirconium (Zr) conversion coating on aluminum using [NEX QC+](#).

## Background

Aluminum and steel are often coated with a protective conversion coating, also called passivate or passivation coating, to prevent oxidation and corrosion of the base metal. Conversion coatings include chromium (Cr), titanium (Ti), vanadium (V), manganese (Mn), nickel (Ni), phosphorus (P), or zirconium (Zr). A phosphate coating may also be applied to minimize wear on cutting tools and stamping machines.


Aluminum is often coated for use in aircraft parts, aluminum window frames, and other similar industries where the aluminum is exposed to weathering. Steel for the automotive industry is typically first galvanized with a zinc coating before the conversion coating is applied. Protected steel is also used for outdoor sheds and other similar uses where steel is exposed to weathering. Conversion coating also helps in the retention of paint for the final finished product.

## Calibration

An empirical calibration was built using a set of standards assayed by careful weigh-strip-weigh.

<b>Element: Zr</b>		
<b>Units: mg/ft<sup>2</sup></b>		
Sample I.D.	Standard value	Calculated value
STD 1	0.98	1.010

STD 2	2.21	2.267
STD 3	3.52	3.446
STD 4	4.81	4.738
STD 5	6.71	6.764
STD 6	15.26	15.265

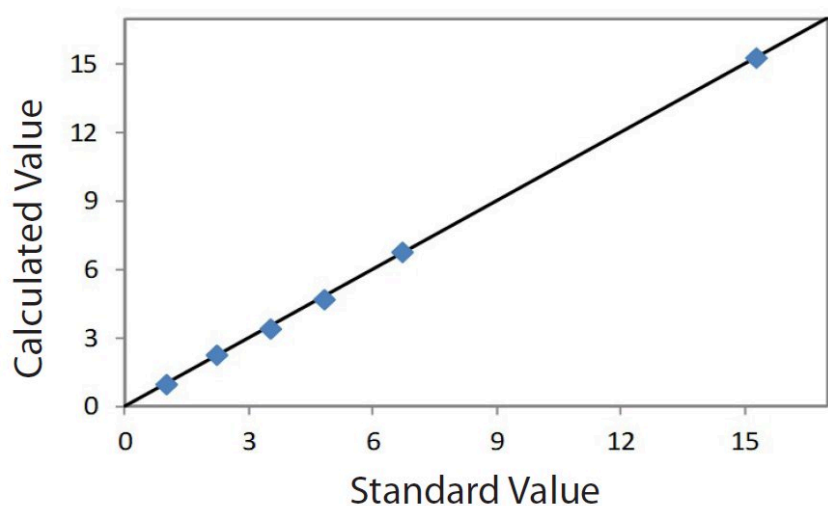

Correlation plot Zr on aluminum

## Recovery and repeatability

To demonstrate repeatability (precision), three calibration standards were chosen. Each sample was measured in a static position for ten repeat analyses with typical results shown below.

<b>Element: Zr</b>				
<b>Units: mg/ft<sup>2</sup></b>				
Sample I.D.	Standard value	Average value	Std. dev	% Relative
STD 1	0.98	1.082	0.029	3.0
STD 4	4.81	4.690	0.037	0.8
STD 6	15.26	15.217	0.024	0.2

## Sample surface and positioning

Aluminum may be surfaced with different patterns to give products various esthetic looks or physical properties. Such surfacing techniques may leave the aluminum smooth, may be brushed aluminum with a directional grain pattern, or may be a hatched or other pattern. The samples demonstrated here had a brushed directional grain pattern.

---

## Conclusion

The performance shown here demonstrates that NEX QC+ provides excellent sensitivity and performance for measuring zirconium conversion coatings on aluminum.

---

## Related products



### NEX QC Series

Combines quality, affordability, and performance for a wide range of applications