

[View on rigaku.com](https://www.rigaku.com)

XRF1117 - Lead analysis in gasoline — ASTM D5059-21 — using WDXRF ZSX Primus IVi

Introduction

Lead (Pb), in the form of alkyl lead such as tetraethyl lead, used to be added to gasoline as an antiknock additive to increase the octane rating. However, alkyl lead is highly poisonous. Gasoline with alkyl lead, called leaded gasoline, causes air pollution. Nowadays, most countries, except in Africa, ban leaded gasoline for vehicles. Alkyl lead is still used as an additive in aviation gasoline.

Although lead-free gasoline, called unleaded gasoline, is common in most countries, lead occurs as a contaminant—either unintentionally or intentionally—in some countries. Therefore, it is necessary to check the lead concentration in gasoline.

This application note demonstrates quantitative analysis of low concentration lead in gasoline according to ASTM D5059-21 on Rigaku ZSX Primus IVi, a wavelength dispersive X-ray fluorescence (WDXRF) spectrometer.

Instrument

The ZSX Primus IVi, a tube-below sequential WDXRF spectrometer, is optimized for the routine analysis that today's petroleum laboratories need to perform. The programmable, switchable vacuum seal between the sample and optical chambers can keep the optical chamber under vacuum with the sample chamber under helium, and, therefore, minimize helium gas consumption and time of atmosphere change in the sample chamber. The spectrometer is equipped with a 4 kW or 3 kW X-ray tube and analyzing crystals covering θ to 2θ in the standard configuration.

The system software is designed for ease of use in routine analyses. The Flowbar in quantitative analysis guides users in establishing calibration. The Sample ID Table and the Program Operation help operators carry out daily analysis.

Measurements were performed on the ZSX Primus IVi with a 4 kW X-ray tube operating at 50 kV and 40 mA using a LiF(200) analyzing crystal and the S2 slit included in the standard configuration. Primary beam filter Al125, inserted between the sample and the X-ray tube, protects the X-ray tube window against damage from falling samples during measurement and also improves peak-to-background ratio.

Analysis method of ASTM D5059-21

The analysis demonstrated in this application note followed Method C of ASTM D5059-21, for low concentration lead, 0.01 g Pb / US gal – 0.5 g Pb / US gal (0.0026 g / L – 0.13 g / L).

In Method C of ASTM D5059-21, the internal standard method with bismuth (Bi) is employed. For each solution of calibration standards or analysis gasoline samples, exactly 20 mL is taken and 2 mL of the Bi internal standard solution (3.00 g Bi / US gal) is added and mixed thoroughly.

X-ray intensities of the following three lines are measured:

- Pb-L α (at 1.175 Å)
- Bi-L α (at 1.144 Å)
- Background (1.194 Å)

The intensity ratio R is determined by the following formula:

$$R = (A - C) / B$$

where A: counting rate at 1.175 Å (Pb-L α)

B: counting rate at 1.144 Å (Bi-L α)

C: counting rate at 1.194 Å (background).

The corrected ratio R_C is determined as follows:

$$R_C = R - R_b$$

where R_b : ratio (R) for blank.

A calibration curve is made with R_C and the slope S of the resulting line is determined as follows:

$$S = (g \text{ Pb} / \text{US gal}) / R_C$$

The lead content of analysis samples is calculated as follows:

$$\text{Pb content (g / US gal)} = S \times R_C$$

The obtained values of the lead content are reported to the nearest 0.005 g / US gal.

Calibration and standards

"Standards for Lead in Gasoline" (isooctane base), which consists of 0.000 (blank), 0.001, 0.005, 0.010, 0.050, 0.100 and 0.300 g Pb / US gal, and "Internal Standards for XRF Analysis; Bi 0.793 g / L" provided by VHG Labs (LGC Standards) were used to make a calibration curve.

For each calibration standard, exactly 20 mL was taken and 2 mL of the Bi internal standard was added and mixed thoroughly. Approximately 6 g of the mixture of each calibration standard and the Bi internal standard was poured into the liquid cell, Chemplex XRF Sample Cups 1095, with 6.0 μm polyester film.

The X-ray intensities of Pb-L α , Bi-L α and the background (1.194 Å) were measured for each of the calibration standards. The counting time was 20 seconds for each line. Then, R_C was calculated and the calibration curve was made. The obtained calibration curve was shown in Figure 1.

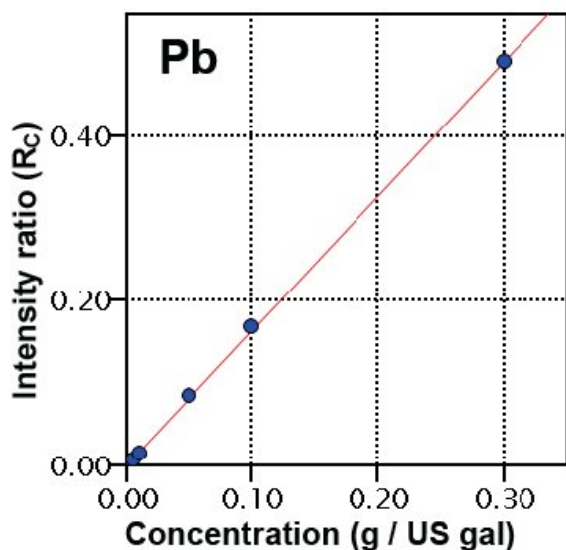


Figure 1: Calibration curve of lead (Pb) in gasoline by bismuth (Bi) internal standard. The accuracy of calibration is 0.0016 g / US gal.

The accuracy of calibration was calculated by the following formula:

$$Accuracy = \sqrt{\frac{\sum_i (C_i - \hat{C}_i)^2}{n-m}}$$

C_i : certified value of standard sample

\hat{C}_i : calculated value of standard sample

n : number of standard samples.

m: degree of freedom (2: linear)

Repeatability test

For repeatability tests, solutions with approximately 0.01 g Pb / US gal and 0.05 g Pb / US gal were prepared by mixing an organometallic standard of Pb with isooctane 75% - toluene 25%.

Repeatability tests were carried out for the above two solutions (~0.01 g Pb / US gal and ~0.05 g Pb / US gal). For the same sample, two aliquots were prepared and quantified with the calibration; this process was repeated twenty times. The test results are tabulated in Table 1, in which the average and the difference of two aliquots each are shown and **r** represents "repeatability" defined in ASTM D5059-21, which states that the difference between successive test results obtained by the same operator with the same apparatus under constant operation conditions on identical test material would, in the long run, in the normal and correct operation of the test method, exceed the following values only in one case in twenty:

$$r = 0.007 + 0.14 X$$

X: g Pb/US gal.

The test results shown in Table 1, where the difference between two aliquots does not exceed the repeatability (**r**), prove that the performance of the ZSX Primus IVI meets the requirement of ASTM D5059-21.

Another test by 10-time measurements was carried out using the same solutions. Ten aliquots were prepared and measured consecutively. The results of the 10-time measurement are tabulated in Table 2.

Table 1: Result of repeatability test for ASTM D5059-21.

Sample: ~0.01 g Pb / US gal				(unit: g Pb / US gal)	
Run	1st aliq.	2nd aliq.	Average	Diff.	r (limit)
01	0.005	0.010	0.008	0.005	0.008
02	0.010	0.010	0.010	0.000	0.008
03	0.010	0.010	0.010	0.000	0.008
04	0.010	0.010	0.010	0.000	0.008
05	0.010	0.005	0.008	0.005	0.008
06	0.005	0.005	0.005	0.000	0.008
07	0.010	0.005	0.008	0.005	0.008
08	0.010	0.010	0.010	0.000	0.008
09	0.010	0.010	0.010	0.000	0.008
10	0.010	0.010	0.010	0.000	0.008
11	0.010	0.010	0.010	0.000	0.008
12	0.010	0.010	0.010	0.000	0.008
13	0.010	0.005	0.008	0.005	0.008
14	0.005	0.010	0.008	0.005	0.008
15	0.010	0.010	0.010	0.000	0.008
16	0.005	0.010	0.008	0.005	0.008
17	0.005	0.010	0.008	0.005	0.008
18	0.010	0.005	0.008	0.005	0.008
19	0.005	0.005	0.005	0.000	0.008
20	0.005	0.010	0.008	0.005	0.008

Sample: ~0.05 g Pb / US gal				(unit: g Pb / US gal)	
Run	1st aliq.	2nd aliq.	Average	Diff.	r (limit)
01	0.040	0.045	0.043	0.005	0.013
02	0.045	0.045	0.045	0.000	0.013
03	0.045	0.040	0.043	0.005	0.013
04	0.040	0.040	0.040	0.000	0.013

05	0.040	0.040	0.040	0.000	0.013
06	0.045	0.040	0.043	0.005	0.013
07	0.045	0.045	0.045	0.000	0.013
08	0.045	0.045	0.045	0.000	0.013
09	0.045	0.045	0.045	0.000	0.013
10	0.045	0.045	0.045	0.000	0.013
11	0.045	0.045	0.045	0.000	0.013
12	0.045	0.045	0.045	0.000	0.013
13	0.045	0.045	0.045	0.000	0.013
14	0.045	0.040	0.043	0.005	0.013
15	0.040	0.040	0.040	0.000	0.013
16	0.040	0.040	0.040	0.000	0.013
17	0.040	0.045	0.043	0.005	0.013
18	0.045	0.045	0.045	0.000	0.013
19	0.045	0.045	0.045	0.000	0.013
20	0.045	0.045	0.045	0.000	0.013

Table 2: Repeatability test results by 10-time measurement.

Sample	~0.01 g Pb / US gal	~0.05 g Pb / US gal
Run	g Pb / US gal	g Pb / US gal
01	0.005	0.040
02	0.005	0.045
03	0.010	0.045
04	0.010	0.045
05	0.010	0.045
06	0.010	0.045
07	0.010	0.045
08	0.010	0.045
09	0.005	0.045
10	0.010	0.040
Average	0.009	0.044
	(0.0024 g / L)	(0.0116 g / L)
Std. Dev.	0.0024	0.0021

	(0.0006 g / L)	(0.0006 g / L)
--	----------------	----------------

Conclusion

Lead in gasoline can be routinely analyzed with high accuracy and precision on the ZSX Primus IVi, a sequential WDXRF spectrometer. This application note demonstrates that the performance of the ZSX Primus IVi meets the requirement of ASTM D5059-21.

Reference

ASTM 5059-21 Standard Test Methods for Lead in Gasoline by X-Ray Spectroscopy

Related products



ZSX Primus IVi

High-power, tube-below, sequential WDXRF spectrometer with new ZSX Guidance expert system software