[rigaku.com](https://www.rigaku.com)で見る

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In order to deconvolute valence density from thermal displacement, experimental charge-density refinements require high-quality data from a vast resolution range. As the atomic scattering factors for high resolution reflections tend to be very small, these reflections are very weak. Therefore, good crystal quality, highly sensitive, noise-free and high-resolution detectors, as well as stable and intense radiation sources are required to make them visible. The low-resolution reflections of such high-quality crystals on the other hand tend to be extremely strong and therefore regularly meet or even exceed the dynamic range of popular phosphor-based integrating detectors. These instrument types require special experimental approaches to overcome the hardware limitations: low resolution data often need to be recorded, by appending additional extremely fast scans or artificially lowering the intensity with attenuators. Using scintillation detectors, the strongest reflections are regularly determined with low redundancy and inferior data quality. In the context of relatively few, low-resolution reflections bearing all information about the anisotropic atomic density, this praxis is sub-optimal. At the same time, scintillation detectors are always affected by electronic noise, ever growing with exposure time and indistinguishable from—and therefore often obscuring—reflection intensities. Both experimental challenges—the determination of extremely weak, as well as extremely strong reflections—can be met by changing from scintillation detectors to direct detection hybrid photon counting detectors. Rigaku implemented this groundbreaking new technology in the HyPix detectors create new possibilities for the collection of high-resolution, high quality data that are required for experimental charge-density.

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The HyPix detector series combines the requirements of experimental charge-density with an excellent spatial resolution of $100\ \mu\text{m}$. As illustrated in Figure 2, very strong and weak reflections are recorded on the same image. The background noise merely stems from stray radiation and air scattering—not from electronic noise as in scintillation detectors. It is therefore zero or close to zero for most parts of the detector (see Figure 2). The HyPix is therefore an excellent choice of detector for experimental charge-density determination.

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