

[View on rigaku.com](https://www.rigaku.com)

BATT1034: Evaluation of crystallite size and grain size distribution of fuel cell materials by in-situ XRD and SAXS

Introduction

In fuel cells, it is known that the average particle size of the Pt catalyst used in the fuel electrode, which is a component material, affects the battery life. Such particle size variation can be evaluated using XRD and SAXS. In this study, the crystallite size of Pt nanoparticles was evaluated by in-situ XRD measurements in a hydrogen gas atmosphere (room temperature) using a corrosion-resistant infrared heated sample high temperature attachment. Small-angle X-ray scattering (SAXS) measurements were also performed before and after the introduction of hydrogen gas to investigate the average particle size and size distribution of Pt nanoparticles.

Particle size and structure analysis

- **Analysis:** Fuel cell
- **Use:** Optimizing electrochemical performance
- **Analyzed materials:** Pt as catalyst
- **Analysis method:** Crystallite size analysis, grain size analysis

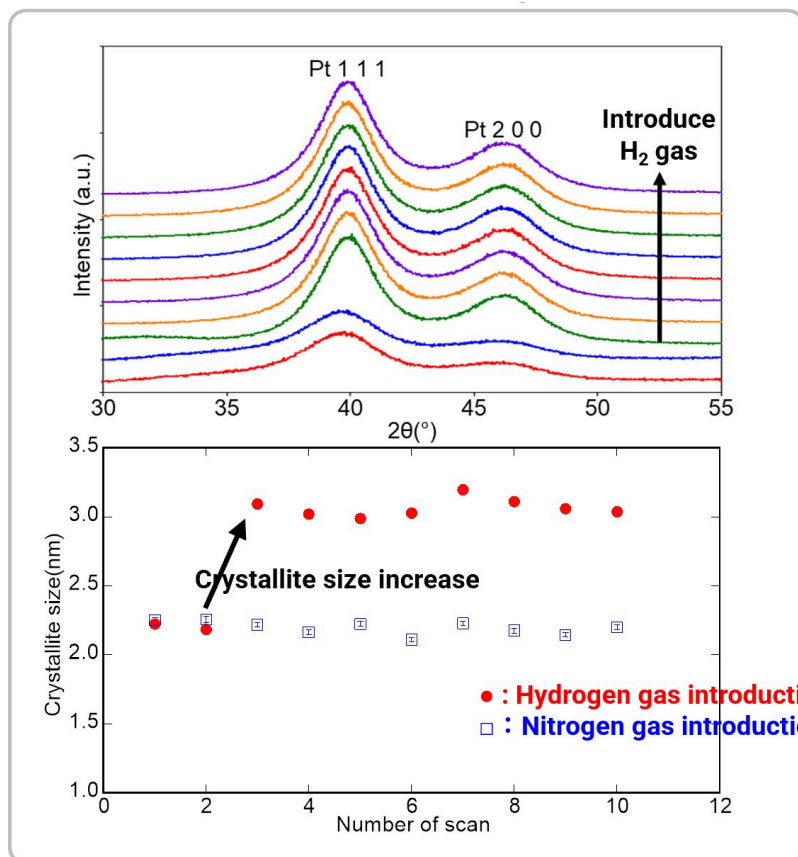


Figure 1: XRD pattern with hydrogen gas introduction and crystallite size change of Pt nanoparticles

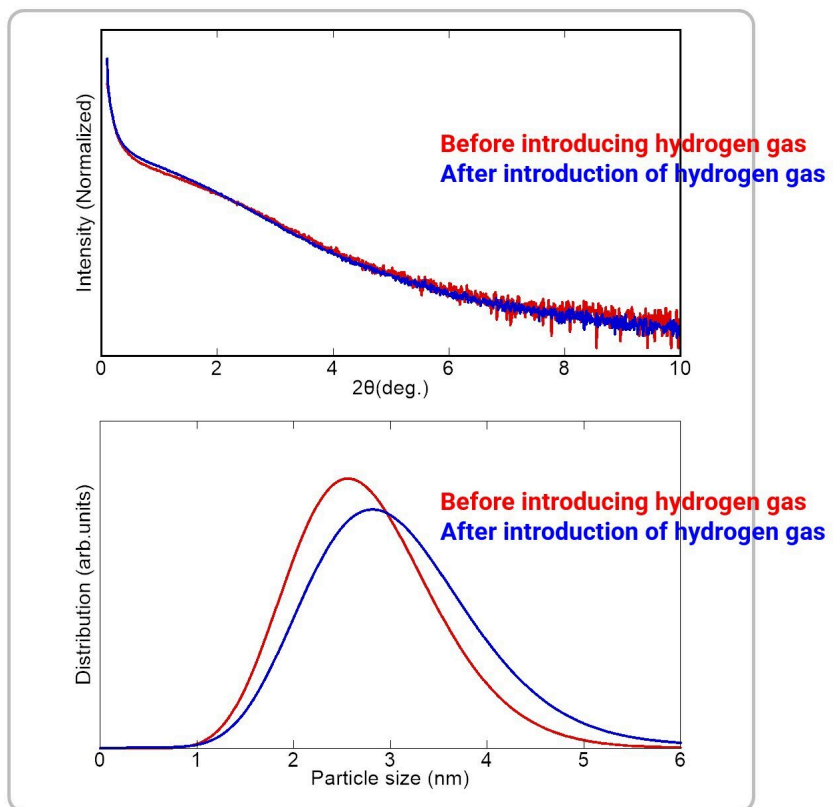


Figure 2: SAXS patterns before and after introduction of hydrogen gas and particle size distribution analysis of Pt nanoparticles

Conclusion

The introduction of hydrogen gas resulted in a sharp change in the diffraction peak shape of the XRD pattern, confirming an increase in the crystallite size of the Pt nanoparticles. The change in the slope of the SAXS profile also suggested a change in particle size. Thus, the ability to evaluate crystallite size and grain size under different environments is a useful method for elucidating the mechanism of structural changes in materials.

Related products



SmartLab

Advanced state-of-the-art high-resolution XRD system powered by Guidance expert system software



SmartLab SE

Highly versatile multipurpose X-ray diffractometer with built-in intelligent guidance



SmartLab Studio II

Windows-based software suite for Rigaku's X-ray diffractometers