

[View on rigaku.com](http://www.rigaku.com)


PHARM019: Quality Confirmation of Disinfectant Solution using 1064 nm Excitation Raman

Introduction

In many cases, manufacturing sites are forced to ship large quantities of hand sanitizers with limited inspection systems due to increasing demand. Cases have been reported in which some of the ingredients do not meet standards or are contaminated with toxic substances such as methanol, which may not be identified by visual or simple inspections. If such risks are overlooked, it may lead to product recalls and loss of trust. With a portable Raman spectrometer, the authenticity of ingredients can be confirmed quickly and nondestructively on site.

Chemical fingerprinting

Analysis:	Pharmaceutical product
Use:	Formulation /Manufacturing (Raw materials)
Analyzed materials:	Hand sanitizer (ethanol)

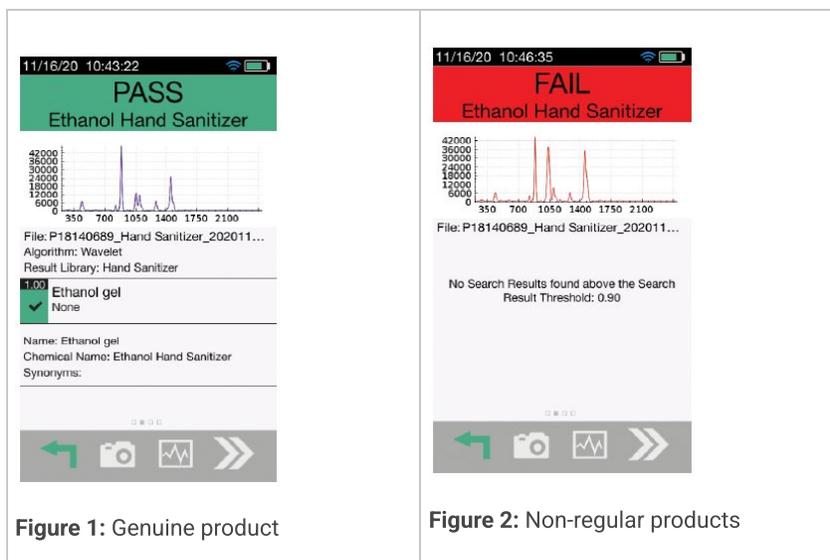

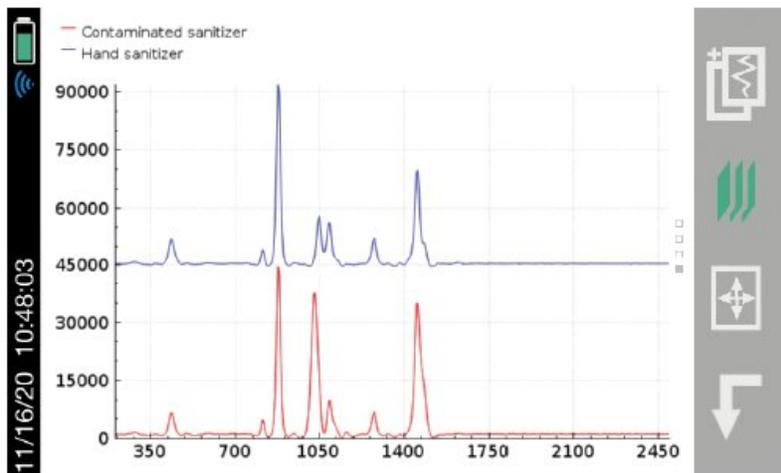


Figure 3: Comparison of Raman spectra
(blue: regular and red: non-regular)

Conclusion

Analysis using a portable Raman spectrometer will show any impurities in a disinfectant. The presence or absence of a substance can be determined immediately on the spot. As shown in Figure 1, the "PASS" mark is displayed for legitimate products, while the "FAIL" mark is displayed for non-legitimate products contaminated with methanol, as shown in Figure 2. As shown in Figure 3, which compares the Raman spectra of the regular product and the non-regular product, a clear peak difference appears around 1000 cm^{-1} . This difference tends to be overlooked by the naked eye. To prevent product recalls and health hazards, a reliable analytical method like this that can be used in the field is indispensable.

Related products



Progeny

Handheld Raman for raw material identification and finished product authentication using 1064 nm Raman analysis.